

In situ HCR v3.0 protocol for samples in solution

This protocol has not been validated for all sample types yet, and should only be used as a template.

Technical support: support@moleculartechnologies.org

Sample pre-treatment

Samples should be prepared in the same manner as for a traditional in situ hybridization, up to the probe hybridization step. This may include permeabilization and protease digestion. Then proceed with the protocol described below.

Multiplexed in situ HCR v3.0 protocol

Detection stage

1. Pre-hybridize samples in 500 μL of 30% probe hybridization buffer for 30 min at 37 °C.
CAUTION: probe hybridization buffer contains formamide, a hazardous material.
2. Prepare probe solution by adding 2 pmol of each probe mixture (odd & even: 1 μL of 2 μM stock per probe mixture) to 500 μL of 30% probe hybridization buffer at 37 °C.
NOTE: For dHCR imaging, use higher concentration of probe to improve probe hybridization efficiency.
3. Remove the pre-hybridization solution and add the probe solution.
4. Incubate samples overnight (12–16 h) at 37 °C.
5. Remove excess probes by washing 4 \times 15 min with 500 μL of 30% probe wash buffer at 37 °C.
CAUTION: probe wash buffer contains formamide, a hazardous material.
NOTE: Wash solutions should be pre-heated to 37 °C before use.
6. Wash samples 3 \times 5 min with 5 \times SSCT at room temperature.

Amplification stage

1. Pre-amplify samples in 500 μL of amplification buffer for 30 min at room temperature.
2. Separately prepare 30 pmol of hairpin H1 and 30 pmol of hairpin H2 by snap cooling 10 μL of 3 μM stock (heat at 95 °C for 90 seconds and cool to room temperature in a dark drawer for 30 min).
NOTE: HCR hairpins H1 and H2 are provided in hairpin storage buffer ready for snap cooling. H1 and H2 should be snap cooled in separate tubes.
3. Prepare hairpin solution by adding all snap-cooled hairpins to 500 μL of amplification buffer at room temperature.
4. Remove the pre-amplification solution and add the hairpin solution.
5. Incubate samples overnight (12–16 h) in the dark at room temperature.
NOTE: For dHCR imaging, amplify for a shorter period of time to ensure single-molecule dots are diffraction-limited.
6. Remove excess hairpins by washing with 500 μL of 5 \times SSCT at room temperature:
 - (a) 2 \times 5 min
 - (b) 2 \times 30 min
 - (c) 1 \times 5 min

Sample mounting for microscopy

Samples can be mounted in SlowFade Gold antifade mountant inside an imaging chamber tailored for the sample of interest. For whole-mount embryos or larvae, a Zeiss 710 NLO inverted confocal microscope equipped with either an LD LCI Plan-Apochromat 25×/0.8 Imm Corr DIC objective or an LD C-Apochromat 40×/1.1 W Korr M27 objective was used in our lab to acquire images.

Buffer recipes for in situ HCR

Probes, amplifiers, probe hybridization buffer, and probe wash buffer should be stored at -20 °C. Amplification buffer should be stored at 4 °C. Keep these reagents on ice at all times during probe and amplifier preparation. Make sure all solutions are well mixed before use.

30% probe hybridization buffer

30% formamide
5× sodium chloride sodium citrate (SSC)
9 mM citric acid (pH 6.0)
0.1% Tween 20
50 µg/mL heparin
1× Denhardt's solution
10% dextran sulfate

For 40 mL of solution

12 mL formamide
10 mL of 20× SSC
360 µL 1 M citric acid, pH 6.0
400 µL of 10% Tween 20
200 µL of 10 mg/mL heparin
800 µL of 50× Denhardt's solution
8 mL of 50% dextran sulfate
Fill up to 40 mL with ultrapure H₂O

30% probe wash buffer

30% formamide
5× sodium chloride sodium citrate (SSC)
9 mM citric acid (pH 6.0)
0.1% Tween 20
50 µg/mL heparin

For 40 mL of solution

12 mL formamide
10 mL of 20× SSC
360 µL 1 M citric acid, pH 6.0
400 µL of 10% Tween 20
200 µL of 10 mg/mL heparin
Fill up to 40 mL with ultrapure H₂O

Amplification buffer

5× sodium chloride sodium citrate (SSC)
0.1% Tween 20
10% dextran sulfate

For 40 mL of solution

10 mL of 20× SSC
400 µL of 10% Tween 20
8 mL of 50% dextran sulfate
Fill up to 40 mL with ultrapure H₂O

5× SSCT

5× sodium chloride sodium citrate (SSC)
0.1% Tween 20

For 40 mL of solution

10 mL of 20× SSC
400 µL of 10% Tween 20
Fill up to 40 mL with ultrapure H₂O

50% dextran sulfate

50% dextran sulfate

For 40 mL of solution

20 g of dextran sulfate powder
Fill up to 40 mL with ultrapure H₂O

Reagents and supplies

Formamide (Deionized) (Ambion Cat. # AM9342)

20× sodium chloride sodium citrate (SSC) (Life Technologies Cat. # 15557-044)

Heparin (Sigma Cat. # H3393)

50% Tween 20 (Life Technologies Cat. # 00-3005)

50× Denhardt's solution (Life Technologies Cat. # 750018)

Dextran sulfate, mol. wt. > 500,000 (Sigma Cat. # D6001)

SlowFade Gold antifade mountant (Life Technologies Cat. # S36936)

Citation Notes

Molecular Technologies (moleculartechnologies.org) is a non-profit academic resource within the Beckman Institute at Caltech. For citation, please select from the list below as appropriate for your application:

- **In situ HCR v3.0**
Automatic background suppression for dramatically enhanced performance (signal-to-background and quantitative precision) and ease-of-use (no probe set optimization for new targets) (Choi *et al.*, 2018). v3.0 supports three quantitative modes: qHCR imaging – analog mRNA relative quantitation with subcellular resolution; qHCR flow cytometry – analog mRNA relative quantitation for high-throughput single-cell analysis; dHCR imaging – digital mRNA absolute quantitation. Protocols for v3.0 in diverse organisms are adapted from the zoo paper.
- **Quantitative HCR (qHCR) imaging**
mRNA relative quantitation with subcellular resolution in the anatomical context of whole-mount vertebrate embryos; read-out/read-in analysis framework (Trivedi *et al.*, 2018).
- **Zoo paper**
Protocols for multiplexed mRNA imaging in diverse sample types (Choi *et al.*, 2016):

bacteria in suspension	whole-mount zebrafish embryos and larvae
whole-mount fruit fly embryos	whole-mount chicken embryos
whole-mount worm larvae	whole-mount mouse embryos
whole-mount sea urchin embryos	FFPE human tissue sections
- **Single-molecule HCR imaging**
Single-molecule mRNA imaging in thick autofluorescent samples (e.g., 0.5 mm adult mouse brain sections) (Shah *et al.*, 2016).
- **Multiplexed quantitative HCR (qHCR) northern blots**
Simultaneous quantification of RNA target size and abundance with signal amplification for up to 5 target RNAs (Schwarzkopf & Pierce, 2016).
- **In situ HCR v2.0**
2nd generation in situ HCR technology (v2.0) using DNA probes and DNA HCR amplifiers: 10× increase in signal, 10× reduction in cost, dramatic increase in reagent durability (Choi *et al.*, 2014).
- **Shielded covalent (SC) probes**
Highly selective covalent capture of DNA and RNA targets, overcoming the longstanding tradeoff between selectivity and durable target capture (Vieregg *et al.*, 2013).
- **In situ HCR v1.0**
1st generation in situ HCR technology (v1.0) using RNA probes and RNA HCR amplifiers: multiplexed mRNA imaging in whole-mount vertebrate embryos with simultaneous signal amplification for up to 5 target mRNAs (Choi *et al.*, 2010).
- **Hybridization chain reaction (HCR) mechanism**
Isothermal enzyme-free molecular signal amplification (Dirks & Pierce, 2004).

We are happy to provide advice and technical support: support@moleculartechnologies.org.

– The Molecular Technologies Team

References

- Choi, H. M. T., Chang, J. Y., Trinh, L. A., Padilla, J. E., Fraser, S. E., & Pierce, N. A. (2010). Programmable in situ amplification for multiplexed imaging of mRNA expression. *Nat Biotechnol*, **28**(11), 1208–12. ([pdf](#), [supp info](#), [supp movie 1](#), [supp movie 2](#))
- Choi, H. M. T., Beck, V. A., & Pierce, N. A. (2014). Next-generation in situ hybridization chain reaction: higher gain, lower cost, greater durability. *ACS Nano*, **8**(5), 4284–4294. ([pdf](#), [supp info](#), [supp movie](#))
- Choi, H. M. T., Calvert, C. R., Husain, N., Huss, D., Barsi, J. C., Deverman, B. E., Hunter, R. C., Kato, M., Lee, S. M., Abelin, A. C. T., Rosenthal, A. Z., Akbari, O. S., Li, Y., Hay, B. A., Sternberg, P. W., Patterson, P. H., Davidson, E. H., Mazmanian, S. K., Prober, D. A., van de Rijn, M., Leadbetter, J. R., Newman, D. K., Readhead, C., Bronner, M. E., Wold, B., Lansford, R., Sauka-Spengler, T., Fraser, S. E., & Pierce, N. A. (2016). Mapping a multiplexed zoo of mRNA expression. *Development*, **143**, 3632–3637. ([pdf](#), [supp info](#), [fruit fly movie](#), [sea urchin movie](#), [zebrafish movie](#), [chicken movie](#), [mouse movie](#))
- Choi, H. M. T., Schwarzkopf, M., Fornace, M. E., Acharya, A., Artavanis, G., Stegmaier, J., Cunha, A., & Pierce, N. A. (2018). Third-generation in situ hybridization chain reaction: multiplexed, quantitative, sensitive, versatile, robust. *Development*, **145**, doi:10.1242/dev.165753. ([pdf](#), [supp info](#), [Dot Analysis 1.0 package](#))
- Dirks, R. M., & Pierce, N. A. (2004). Triggered amplification by hybridization chain reaction. *Proc Natl Acad Sci USA*, **101**(43), 15275–15278. ([pdf](#))
- Schwarzkopf, M., & Pierce, N. A. (2016). Multiplexed miRNA northern blots via hybridization chain reaction. *Nucleic Acids Res*, **44**(15), e129. ([pdf](#), [supp info](#))
- Shah, S., Lubeck, E., Schwarzkopf, M., He, T.-F., Greenbaum, A., Sohn, C. H., Lignell, A., Choi, H. M. T., Gradi-naru, V., Pierce, N. A., & Cai, L. (2016). Single-molecule RNA detection at depth via hybridization chain reaction and tissue hydrogel embedding and clearing. *Development*, **143**, 2862–2867. ([pdf](#), [supp info](#), [supp movie 1](#), [supp movie 2](#), [supp movie 3](#))
- Trivedi, V., Choi, H. M. T., Fraser, S. E., & Pierce, N. A. (2018). Multidimensional quantitative analysis of mRNA expression within intact vertebrate embryos. *Development*, **145**, doi:10.1242/dev.156869. ([pdf](#), [supp info](#), [Read-out/Read-in 1.0 package](#))
- Vieregg, J. R., Nelson, H. M., Stoltz, B. M., & Pierce, N. A. (2013). Selective nucleic acid capture with shielded covalent probes. *J Am Chem Soc*, **135**(26), 9691–9699. ([pdf](#), [supp info](#))