

## In situ HCR v3.0 protocol for whole-mount zebrafish larvae (*Danio rerio*)

This protocol has been optimized for embryos at 27 hpf and larvae at 5 dpf. Other developmental stages may require additional optimization.

Technical support: [support@moleculartechnologies.org](mailto:support@moleculartechnologies.org)

### Sample preparation

1. Collect zebrafish embryos and incubate at 28 °C in a petri dish with egg H<sub>2</sub>O.
2. Exchange egg H<sub>2</sub>O with egg H<sub>2</sub>O containing 0.003% of 1-phenyl 2-thiourea (PTU) when embryos reach 12 hpf.  
*NOTE: PTU inhibits melanogenesis but can be toxic at high concentrations. PTU treatment must start before the initial pigmentation occurs as PTU does not remove pigment that has already formed. PTU treatment is not necessary for 27 hpf (hours post-fertilization) embryos.*
3. Replace with fresh egg H<sub>2</sub>O containing 0.003% of PTU everyday until the larvae reach 5 dpf (days post-fertilization).
4. Transfer ~40 larvae (5 dpf) to a 2 mL eppendorf tube and remove excess egg H<sub>2</sub>O.
5. Fix larvae in 2 mL of 4% paraformaldehyde (PFA) for 24 h at 4 °C.  
*CAUTION: Use PFA with extreme care as it is a hazardous material.*  
*NOTE: Use fresh PFA and cool to 4 °C before use to avoid increased autofluorescence.*
6. Wash larvae 3 × 5 min with 1 mL of 1× phosphate-buffered saline (PBS) to stop the fixation.  
*NOTE: Avoid using calcium chloride and magnesium chloride in PBS as this leads to increased autofluorescence in the samples.*
7. Dehydrate and permeabilize with a series of methanol (MeOH) washes (1 mL each):
  - (a) 100% MeOH for 4 × 10 min
  - (b) 100% MeOH for 1 × 50 min.
8. Store larvae at -20 °C overnight before use.  
*NOTE: Larvae can be stored for six months at -20 °C.*

## Buffer recipes for sample preparation

### 6% PTU stock solution

6% PTU

### For 100 mL of solution

6 g of 1-phenyl 2-thiourea powder

Fill up to 100 mL with egg H<sub>2</sub>O

Heat solution at 50–60 °C overnight to dissolve powder

### 0.3% PTU in egg H<sub>2</sub>O

0.3% PTU

### For 50 mL of solution

2.5 mL of 6% PTU

Fill up to 50 mL with with egg H<sub>2</sub>O

### 4% Paraformaldehyde (PFA)

4% PFA

1× PBS

### For 25 mL of solution

1 g of PFA powder

25 mL of 1× PBS

Heat solution at 50–60 °C to dissolve powder

*NOTE: Avoid using calcium chloride and magnesium chloride in PBS as this leads to increased autofluorescence in the samples.*

## Multiplexed in situ HCR v3.0 protocol

### Detection stage

1. Transfer the required number of larvae for an experiment to a 2 mL eppendorf tube.
2. Rehydrate with a series of graded 1 mL MeOH/PBST washes for 5 min each at room temperature:
  - (a) 75% MeOH / 25% PBST
  - (b) 50% MeOH / 50% PBST
  - (c) 25% MeOH / 75% PBST
  - (d) 5 × 100% PBST.
3. Treat 5 dpf larvae with 1 mL of proteinase K (30 µg/mL) for 45 min at room temperature.  
*NOTE: Proteinase K concentration and treatment time should be reoptimized for each batch of proteinase K, or for samples at a different developmental stage. Skip proteinase K treatment and postfixation (steps 3–6) for embryos 30 hpf and younger.*
4. Wash larvae two times with PBST (1 mL each) without incubation.
5. Postfix with 1 mL of 4 % PFA for 20 min at room temperature.
6. Wash larvae 5 × 5 min with 1 mL of PBST.
7. For each sample, move 8 larvae to a 1.5 mL eppendorf tube.
8. Pre-hybridize with 350 µL of 30% probe hybridization buffer for 30 min at 37 °C.  
*CAUTION: probe hybridization buffer contains formamide, a hazardous material.*
9. Prepare probe solution by adding 2 pmol of each probe mixture (odd & even: 1 µL of 2 µM stock per probe mixture) to 500 µL of probe hybridization buffer at 37 °C.  
*NOTE: For dHCR imaging use higher probe concentration to improve probe hybridization efficiency.*
10. Remove the pre-hybridization solution and add the probe solution.
11. Incubate larvae overnight (12–16 h) at 37 °C.
12. Remove excess probes by washing embryos 4 × 15 min with 500 µL of 30% probe wash buffer at 37 °C.  
*CAUTION: probe wash buffer contains formamide, a hazardous material.*  
*NOTE: Wash solutions should be pre-heated to 37 °C before use.*
13. Wash embryos 3 × 5 min with 5× SSCT at room temperature.

### Amplification stage

1. Pre-amplify larvae with 350  $\mu\text{L}$  of amplification buffer for 30 min at room temperature.
2. Separately prepare 30 pmol of hairpin H1 and 30 pmol of hairpin H2 by snap cooling 10  $\mu\text{L}$  of 3  $\mu\text{M}$  stock (heat at 95  $^{\circ}\text{C}$  for 90 seconds and cool to room temperature in a dark drawer for 30 min).  
*NOTE: HCR hairpins H1 and H2 are provided in hairpin storage buffer ready for snap cooling. H1 and H2 should be snap cooled in separate tubes.*
3. Prepare hairpin solution by adding all snap-cooled hairpins to 500  $\mu\text{L}$  of amplification buffer at room temperature.
4. Remove the pre-amplification solution and add the hairpin solution.
5. Incubate larvae overnight (12–16 h) in the dark at room temperature.  
*NOTE: For dHCR imaging, amplify for a shorter period of time to ensure single-molecule dots are diffraction-limited.*
6. Remove excess hairpins by washing with 500  $\mu\text{L}$  of 5 $\times$  SSCT at room temperature:
  - (a) 2  $\times$  5 min
  - (b) 2  $\times$  30 min
  - (c) 1  $\times$  5 min.

### Sample mounting for microscopy

1. A chamber for mounting the larvae was made by aligning two stacks of Scotch tape (8 pieces per stack) 1 cm apart on a 25 mm  $\times$  75 mm glass slide.
2. Approximately 200  $\mu\text{L}$  of 3% methyl cellulose mounting medium was added between the tape stacks on the slide.
3. Place larvae on the medium and orient for dorsal imaging.
4. Place a 22 mm  $\times$  22 mm No. 1 coverslip on top of the stacks to close the chamber.
5. A Zeiss 710 NLO inverted confocal microscope equipped with an LD C-Apochromat 40 $\times$ /1.1 W Korr M27 objective was used in our lab to acquire zebrafish images.

## Buffer recipes for in situ HCR v3.0

Probes, amplifiers, probe hybridization buffer, and probe wash buffer should be stored at -20 °C. Amplification buffer should be stored at 4 °C. Keep these reagents on ice at all times during probe and amplifier preparation. Make sure all solutions are well mixed before use.

### PBST

1× PBS  
0.1% Tween 20

### For 50 mL of solution

5 mL of 10× PBS  
500 μL of 10% Tween 20  
Fill up to 50 mL with ultrapure H<sub>2</sub>O

### Proteinase K solution

30 μg/mL proteinase K

### For 1 mL of solution

1.5 μL of 20 mg/mL proteinase K  
Fill up to 1 mL with PBST

### Probe hybridization buffer

30% formamide  
5× sodium chloride sodium citrate (SSC)  
9 mM citric acid (pH 6.0)  
0.1% Tween 20  
50 μg/mL heparin  
1× Denhardt's solution  
10% dextran sulfate

### For 40 mL of solution

12 mL formamide  
10 mL of 20× SSC  
360 μL 1 M citric acid, pH 6.0  
400 μL of 10% Tween 20  
200 μL of 10 mg/mL heparin  
800 μL of 50× Denhardt's solution  
8 mL of 50% dextran sulfate  
Fill up to 40 mL with ultrapure H<sub>2</sub>O

### 30% Probe wash buffer

30% formamide  
5× sodium chloride sodium citrate (SSC)  
9 mM citric acid (pH 6.0)  
0.1% Tween 20  
50 μg/mL heparin

### For 40 mL of solution

12 mL formamide  
10 mL of 20× SSC  
360 μL 1 M citric acid, pH 6.0  
400 μL of 10% Tween 20  
200 μL of 10 mg/mL heparin  
Fill up to 40 mL with ultrapure H<sub>2</sub>O

### Amplification buffer

5× sodium chloride sodium citrate (SSC)  
0.1% Tween 20  
10% dextran sulfate

### For 40 mL of solution

10 mL of 20× SSC  
400 μL of 10% Tween 20  
8 mL of 50% dextran sulfate  
Fill up to 40 mL with ultrapure H<sub>2</sub>O

### 5× SSCT

5× sodium chloride sodium citrate (SSC)  
0.1% Tween 20

### For 40 mL of solution

10 mL of 20× SSC  
400 μL of 10% Tween 20  
Fill up to 40 mL with ultrapure H<sub>2</sub>O

### 50% dextran sulfate

50% dextran sulfate

### For 40 mL of solution

20 g of dextran sulfate powder  
Fill up to 40 mL with ultrapure H<sub>2</sub>O

### 3% methyl cellulose

3% methyl cellulose

### For 100 mL of solution

3 g of methyl cellulose powder  
Fill up to 100 mL with ultrapure H<sub>2</sub>O  
Stir overnight at 4 °C with a stir bar to dissolve powder

## **Reagents and supplies**

Paraformaldehyde (PFA) (Sigma Cat. # P6148)  
Methanol (Mallinckrodt Chemicals Cat. # 3016-16)  
Proteinase K, molecular biology grade (NEB Cat. # P8107S)  
Formamide (Deionized) (Ambion Cat. # AM9342)  
20× sodium chloride sodium citrate (SSC) (Life Technologies Cat. # 15557-044)  
50% Tween 20 (Life Technologies Cat. # 00-3005)  
Heparin (Sigma Cat. # H3393)  
50× Denhardt's solution (Life Technologies Cat. # 750018)  
Dextran sulfate, mol. wt. > 500,000 (Sigma Cat. # D6001)  
25 mm × 75 mm glass slide (VWR Cat. # 48300-025)  
22 mm × 22 mm No. 1 coverslip (VWR Cat. # 48366-067)  
Methyl cellulose (Sigma Cat. # M0387)

## Citation Notes

Molecular Technologies ([moleculartechnologies.org](http://moleculartechnologies.org)) is a non-profit academic resource within the Beckman Institute at Caltech. For citation, please select from the list below as appropriate for your application:

- **In situ HCR v3.0**  
Automatic background suppression for dramatically enhanced performance (signal-to-background and quantitative precision) and ease-of-use (no probe set optimization for new targets) (Choi *et al.*, 2018). v3.0 supports three quantitative modes: qHCR imaging – analog mRNA relative quantitation with subcellular resolution; qHCR flow cytometry – analog mRNA relative quantitation for high-throughput single-cell analysis; dHCR imaging – digital mRNA absolute quantitation. Protocols for v3.0 in diverse organisms are adapted from the zoo paper.
- **Quantitative HCR (qHCR) imaging**  
mRNA relative quantitation with subcellular resolution in the anatomical context of whole-mount vertebrate embryos; read-out/read-in analysis framework (Trivedi *et al.*, 2018).
- **Zoo paper**  
Protocols for multiplexed mRNA imaging in diverse sample types (Choi *et al.*, 2016):
 

bacteria in suspension	whole-mount zebrafish embryos and larvae
whole-mount fruit fly embryos	whole-mount chicken embryos
whole-mount worm larvae	whole-mount mouse embryos
whole-mount sea urchin embryos	FFPE human tissue sections
- **Single-molecule HCR imaging**  
Single-molecule mRNA imaging in thick autofluorescent samples (e.g., 0.5 mm adult mouse brain sections) (Shah *et al.*, 2016).
- **Multiplexed quantitative HCR (qHCR) northern blots**  
Simultaneous quantification of RNA target size and abundance with signal amplification for up to 5 target RNAs (Schwarzkopf & Pierce, 2016).
- **In situ HCR v2.0**  
2nd generation in situ HCR technology (v2.0) using DNA probes and DNA HCR amplifiers: 10× increase in signal, 10× reduction in cost, dramatic increase in reagent durability (Choi *et al.*, 2014).
- **Shielded covalent (SC) probes**  
Highly selective covalent capture of DNA and RNA targets, overcoming the longstanding tradeoff between selectivity and durable target capture (Vieregg *et al.*, 2013).
- **In situ HCR v1.0**  
1st generation in situ HCR technology (v1.0) using RNA probes and RNA HCR amplifiers: multiplexed mRNA imaging in whole-mount vertebrate embryos with simultaneous signal amplification for up to 5 target mRNAs (Choi *et al.*, 2010).
- **Hybridization chain reaction (HCR) mechanism**  
Isothermal enzyme-free molecular signal amplification (Dirks & Pierce, 2004).

We are happy to provide advice and technical support: [support@moleculartechnologies.org](mailto:support@moleculartechnologies.org).

– The Molecular Technologies Team

## References

- Choi, H. M. T., Chang, J. Y., Trinh, L. A., Padilla, J. E., Fraser, S. E., & Pierce, N. A. (2010). Programmable in situ amplification for multiplexed imaging of mRNA expression. *Nat Biotechnol*, **28**(11), 1208–12. ([pdf](#), [supp info](#), [supp movie 1](#), [supp movie 2](#))
- Choi, H. M. T., Beck, V. A., & Pierce, N. A. (2014). Next-generation in situ hybridization chain reaction: higher gain, lower cost, greater durability. *ACS Nano*, **8**(5), 4284–4294. ([pdf](#), [supp info](#), [supp movie](#))
- Choi, H. M. T., Calvert, C. R., Husain, N., Huss, D., Barsi, J. C., Deverman, B. E., Hunter, R. C., Kato, M., Lee, S. M., Abelin, A. C. T., Rosenthal, A. Z., Akbari, O. S., Li, Y., Hay, B. A., Sternberg, P. W., Patterson, P. H., Davidson, E. H., Mazmanian, S. K., Prober, D. A., van de Rijn, M., Leadbetter, J. R., Newman, D. K., Readhead, C., Bronner, M. E., Wold, B., Lansford, R., Sauka-Spengler, T., Fraser, S. E., & Pierce, N. A. (2016). Mapping a multiplexed zoo of mRNA expression. *Development*, **143**, 3632–3637. ([pdf](#), [supp info](#), [fruit fly movie](#), [sea urchin movie](#), [zebrafish movie](#), [chicken movie](#), [mouse movie](#))
- Choi, H. M. T., Schwarzkopf, M., Fornace, M. E., Acharya, A., Artavanis, G., Stegmaier, J., Cunha, A., & Pierce, N. A. (2018). Third-generation in situ hybridization chain reaction: multiplexed, quantitative, sensitive, versatile, robust. *Development*, **145**, doi:10.1242/dev.165753. ([pdf](#), [supp info](#), [Dot Analysis 1.0 package](#))
- Dirks, R. M., & Pierce, N. A. (2004). Triggered amplification by hybridization chain reaction. *Proc Natl Acad Sci USA*, **101**(43), 15275–15278. ([pdf](#))
- Schwarzkopf, M., & Pierce, N. A. (2016). Multiplexed miRNA northern blots via hybridization chain reaction. *Nucleic Acids Res*, **44**(15), e129. ([pdf](#), [supp info](#))
- Shah, S., Lubeck, E., Schwarzkopf, M., He, T.-F., Greenbaum, A., Sohn, C. H., Lignell, A., Choi, H. M. T., Gradi-naru, V., Pierce, N. A., & Cai, L. (2016). Single-molecule RNA detection at depth via hybridization chain reaction and tissue hydrogel embedding and clearing. *Development*, **143**, 2862–2867. ([pdf](#), [supp info](#), [supp movie 1](#), [supp movie 2](#), [supp movie 3](#))
- Trivedi, V., Choi, H. M. T., Fraser, S. E., & Pierce, N. A. (2018). Multidimensional quantitative analysis of mRNA expression within intact vertebrate embryos. *Development*, **145**, doi:10.1242/dev.156869. ([pdf](#), [supp info](#), [Read-out/Read-in 1.0 package](#))
- Vieregg, J. R., Nelson, H. M., Stoltz, B. M., & Pierce, N. A. (2013). Selective nucleic acid capture with shielded covalent probes. *J Am Chem Soc*, **135**(26), 9691–9699. ([pdf](#), [supp info](#))